


FLARE Wide Pore C18 Column: Separation of Antibody Drug Conjugate (ADC) Digested Sample

Brentuximab-Vedotin was digested with endoproteinase Lys-C

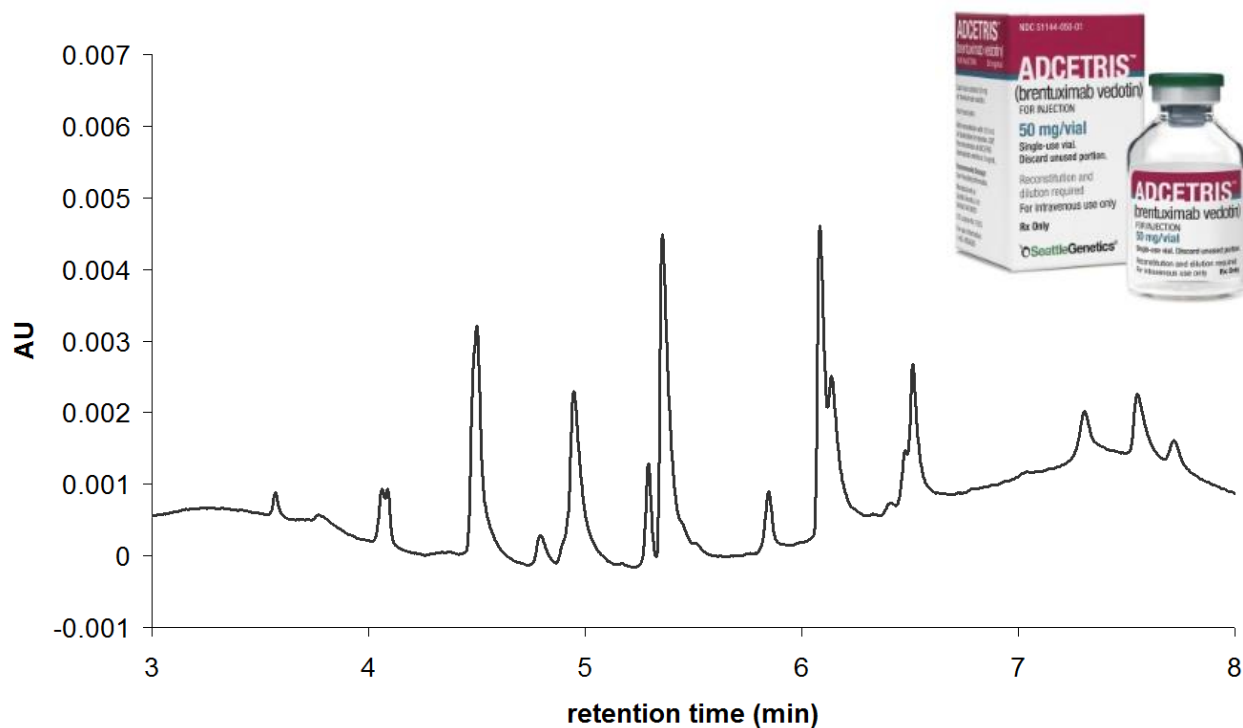

Column Name: FLARE WP C18

Column Dimensions: 100 x 2.1mm (ID 16636.1-2)

Mobile Phase: A: 0.5% TFA; B: 0.5% TFA in ACN

Gradient: 0 – 50% B in 12 mins

Flow Rate: 0.2 mL/min

Temperature: 60 °C

Detection: UV @ 280nm, 20 Hz sampling rate